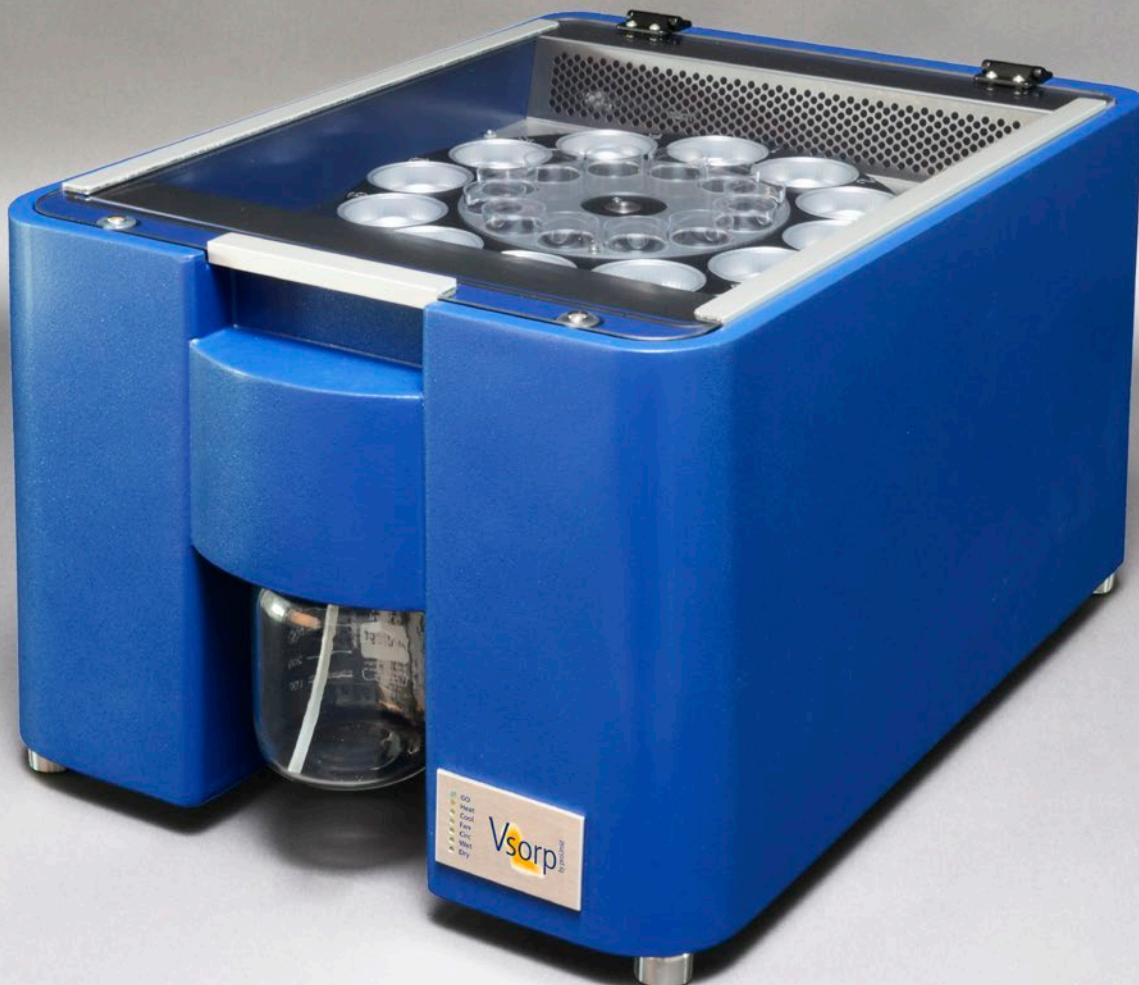
# Dynamic Water Vapor Sorption

proUmid<sup>®</sup>  
GmbH & Co. KG

## Multisample Moisture Sorption Analyzer – Vsorp

# The smart sorption solution

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### High performance

The Vsorp is an easy to operate, high throughput instrument for routine measurements in research, development and quality assurance.

### Advanced design

The Vsorp stands for the combination of proven and reliable components with an advanced design for high reliability, minimized instrument care and low operating cost.

### Value for money

Combining our longstanding experience in the construction of measurement devices with state-of-the-art manufacturing, the Vsorp covers the requirements of an advanced multi-sample instrument at an attractive price.

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# Working with the Vsorp instruments

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### Sample preparation

Sample material is applied directly into the pans on the sample carousel. The easily accessible measurement chamber enables a convenient sample preparation.

### Fully automated process

Measurement and data acquisition are fully automated. The large transparent cover enables the observation of the samples. Acquired data can be exported during a running measurement.

### Operating software

- User-friendly and intuitive interface
- Programmable sorption/desorption cycles and unlimited number of cycle repetitions
- Humidity cycles as step profiles, ramps or plateaus
- Templates for standard measurements
- Optional 21 CFR part 11 compliant software package

### Data export and further evaluation

- Import into an Excel® template with fully configurable graphical and tabulated data
- LIMS compatible data format
- Optional software package for data import and advanced data analysis

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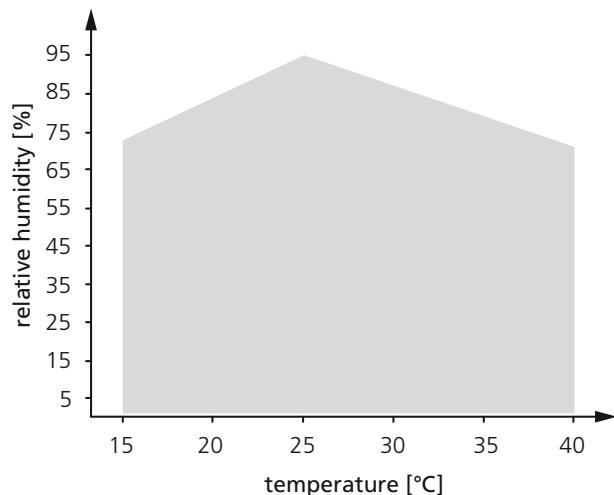
Vsorp instrument	Vsorp Enhanced	Vsorp Plus	Vsorp Basic
Number of samples & sample size	23 samples Ø 33 mm 11 samples Ø 50 mm	11 samples Ø 50 mm	5 samples Ø 86 mm
Temperature range	15 °C to 40 °C ( $\pm 0.1$ K)		
Humidity range*	0 % RH to 95 % RH ( $\pm 0.5$ % RH at 10...30 °C)		
Load range min to max	<30 mg to 220 g	<50 mg to 220 g	<500 mg to 220 g
Balance resolution	1 µg (up to 22 g) 10 µg (up to 220 g)	10 µg (up to 111 g) 100 µg (up to 220 g)	100 µg
Repeatability**	±5 µg (up to 22 g) ±20 µg (up to 220 g)	±20 µg (up to 111 g) ±100 µg (up to 220 g)	±100 µg
Hardware options	Water vapor permeability rate Large objects sample tray		
Software options	21 CFR Part 11 compliant software package Analysis software		

\* Max RH value. See diagram for possible T / RH combinations.

\*\* Root Mean Square (RMS). The specified values assume that the system is installed in an environment suitable for the operation of micro-balances.

Humidity and temperature range for possible T / RH combinations  
at 22 °C ambient temperature.

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# Vsorp special features

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### Humidity & temperature range

The Vsorp is designed for optimized temperature and humidity stability and best performance in the extended room temperature range.

### Proven load cell technology

Baseline stability, accuracy and reproducibility are on a level beyond comparable single sample instruments.

### Exchangeable sample trays

Different sample trays with dish size range from Ø 33–86 mm, as well as pans for large rectangular samples enable the analysis of powders and granules as well as larger items such as packaging or concrete.

### Permeability testing

The optional Permeability Kit enables the investigation of the water vapor transmission rate (WVTR) through films, foils and packaging materials.

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### ProUmid GmbH & Co. KG